

Handling Instruction for Dry Shipper (LN2)

Inserting the Product (For Shipper)

1. Remove the Outer Lid (1) of the Outer Pack
2. Open the Inner Lid (2) of the Tank
3. Take out the Rack (3) from inside the Tank
4. Insert your cryo box(es) into the Rack
5. Insert the Rack into the Tank
6. Close the Inner Lid
7. **LOGGER CHECK:**
 - Please check if the Logger is started
 - If Logger is not started:
 - **PLEASE START LOGGER BY PRESSING AND HOLDING THE START BUTTOM FOR 2 SEC**




Logger is started!



Logger is NOT started!



Press and hold for 2 seconds to start the logger

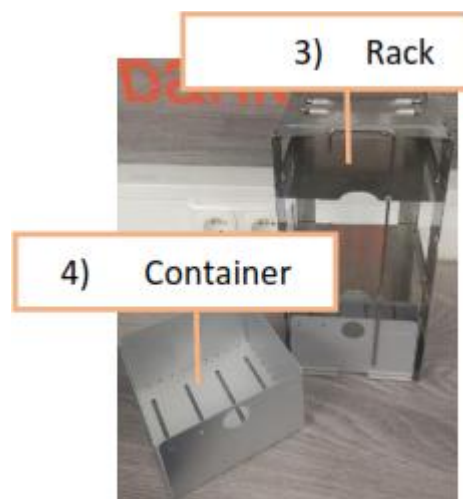

Handling Instruction for Dry Shipper (LN2)

Removing the Product (For Receiver)

1. Remove the Outer Lid (1) of the Outer Pack
2. Open the Inner Lid (2) of the Tank
3. Take out the Rack (3) from inside the Tank
4. Remove your cryo box(es) from the Rack
5. Insert the Rack into the Tank
6. Close the Inner Lid
7. Close the Outer Lid

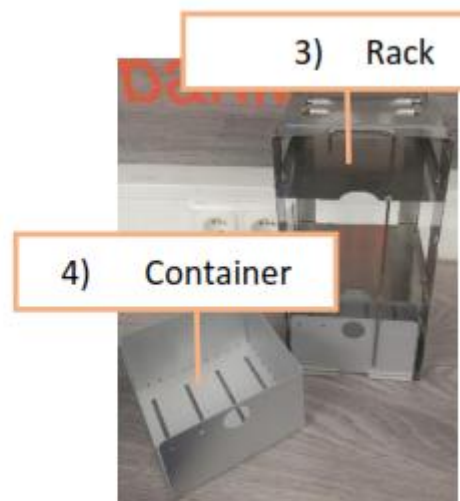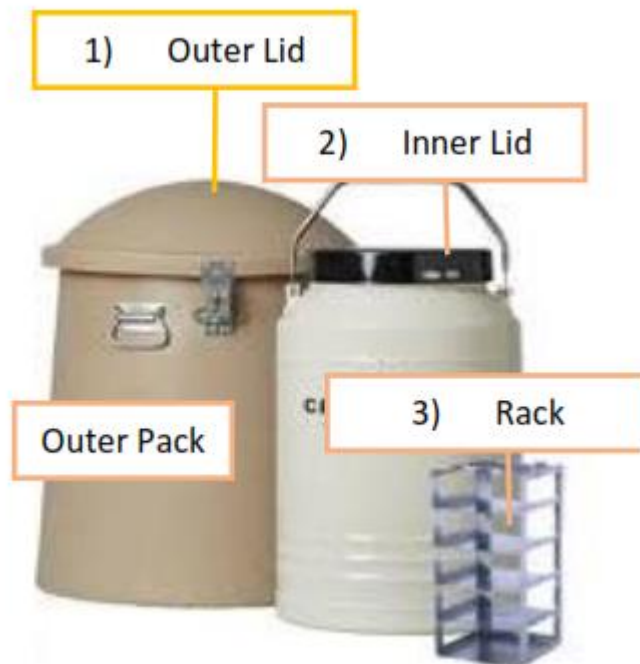


Handling Instruction for Dry Shipper (LN2)

Temperature Logger (For Receivers only)

1. Remove the Logger (1) from the Logger Holder (3)
2. Remove the Temp. Sensor (2) by pulling from the Logger
3. Connect Logger via USB (4) to any PC or Laptop
4. After ~30-40 seconds a folder appears on the desktop with a PDF file inside
5. Save the PDF file
6. Remove Logger from PC/Laptop
7. Connect the Temp. Sensor by pushing to the Logger
8. Insert Logger into Logger Holder



1) Logger

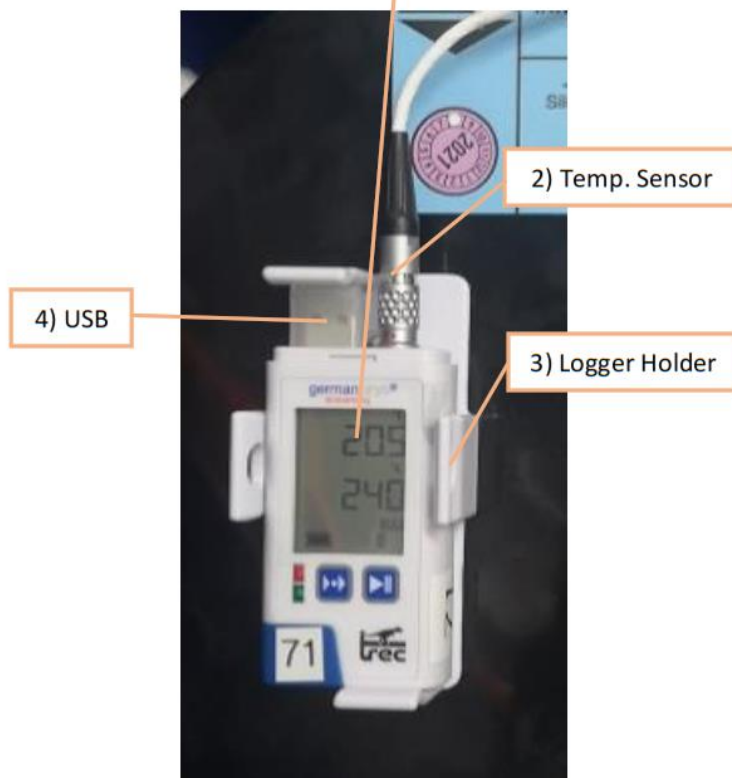

2) Temp. Sensor

4) USB

3) Logger Holder