

Leukapheresis Shipment Instructions

Your facility has been equipped with a specific Thermal box for the transport of Leukapheresis samples to immatics biotechnologies GmbH.

Please adhere to following instructions for sample shipping:

PRE-CONDITIONING OF THERMAL BOX

In order to ensure correct transportation temperature, the thermal box needs to be pre-conditioned before every shipment:

1. Open the box and open the top flap of the cooling frames.
2. Keep the temperature logger and the temperature elements inside.
3. **Store the opened box at 20°C-25°C for a minimum of 24hrs**



PREPARE FOR PICKUP

1. Place Leukapheresis bag(s) into the container
2. Take out the temperature logger and start it by pressing the "Transit" button for at least 3 seconds.
3. Once started, the logger display will show "Transit" on the Screen
4. Place the started logger next to the sample bag(s)
5. Close the transport box and hand it over to the courier

